Quanti-Tray*/2000

Insert and Most Probable Number (MPN) Table

IDX 33/02 – 06/12
WATER ANALYSIS METHODS
www.afnor-validation.org

Quanti-Tray* Certificate of Sterility

This certifies that the enclosed Quanti-Tray/2000 trays have been sterilized with ethylene oxide.

For technical support, please call:

North/South America: +1 207 556 4496 or 1 800 321 0207

Europe: +00800 4339 9111

UK: +44 01638 676800

China: +86 21 61279528

Japan: +81 422 71 5921

Australia: +1800 443 399

idexx.com/water

The method Colilert*-18/Quanti-Tray*/2000 for water analysis is granted NF Validation by AFNOR Certification as an alternative method to the standard ISO 9308-3 for detection and enumeration of *Escherichia coli* β -glucuronidase positive in bathing water, under the Certificate number: IDX 33/02 – 06/12 For more information about end of validity, please refer to the certificate NF Validation available on website mentioned above

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*Quanti-Tray and Colilert are trademarks or registered trademarks of IDEXX Laboratories, Inc. or its affiliates in the United States and/or other countries.

Patent information: idexx.com/patents.

IDEXX

One IDEXX Drive
Westbrook, Maine 04092 USA

Quanti-Tray*/2000

English

Introduction


IDEXX Quanti-Tray*/2000 is designed to give quantitated bacterial counts of 100 mL samples using IDEXX reagent products. Add the reagent/sample mixture to a Quanti-Tray/2000, seal it in a Quanti-Tray* Sealer and incubate per the reagent instructions. Count the number of positive large and small wells and use the Most Probable Number (MPN) Table attached to determine the MPN.

Contents


This package contains sterile Quanti-Tray/2000 trays.

User Instructions


1. Use one hand to hold a Quanti-Tray* upright with the well side facing the palm.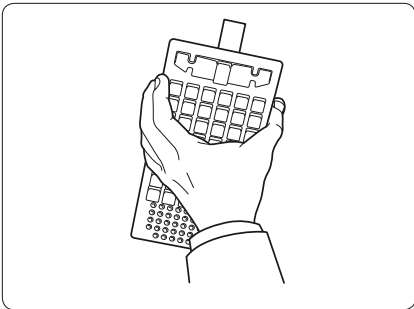


2. Squeeze the upper part of the Quanti-Tray so that the Quanti-Tray bends toward the palm.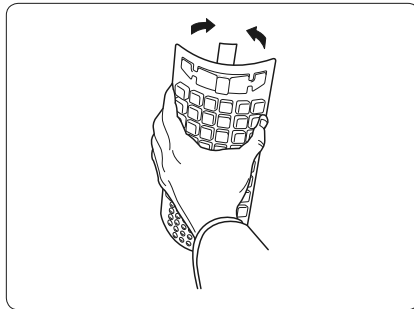


3. Gently pull foil tab to separate the foil from the tray. Avoid touching the inside of the foil or tray.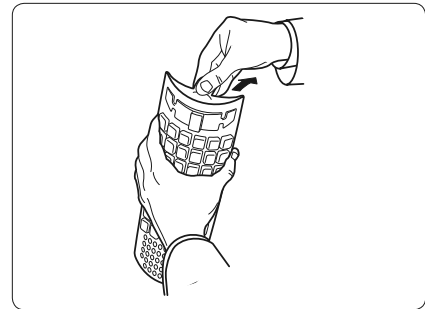

4. Pour the reagent/sample mixture directly into the Quanti-Tray, avoiding contact with the foil tab. Tap the small wells 2–3 times to release any air bubbles. Allow foam to settle.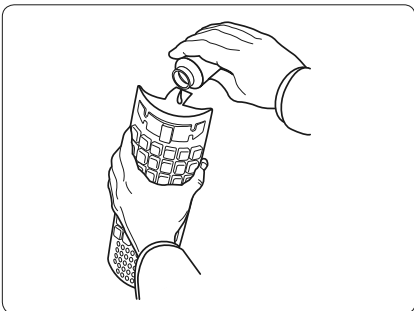

5. Place the sample-filled Quanti-Tray onto the Quanti-Tray/2000 rubber insert of the Quanti-Tray Sealer with the well side (plastic) of the Quanti-Tray facing down.

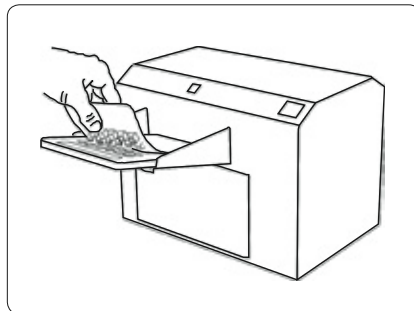6. Seal according to the Quanti-Tray Sealer instructions.
7. Incubate according to reagent instructions.
8. Count large and small positive wells and refer to the Quanti-Tray/2000 MPN table to find the MPN.[†]
9. Dispose of media in accordance with good laboratory practices.

For technical support, please call:

North/South America: +1 207 556 4496 or 1 800 321 0207

Europe: +00800 4339 9111

UK: +44 01638 676800

China: +86 21 61279528

Japan: +81 422 71 5921

Australia: +1800 443 399

idexx.com/water

[†]Download the IDEXX MPN generator software for automated Quanti-Tray results at idexx.com/mpngenerator.

*Quanti-Tray is a trademark or registered trademark of IDEXX Laboratories, Inc. in the United States and/or other countries.

© 2013 IDEXX Laboratories, Inc. All rights reserved.

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